# ACOMP: The Automated system that allows continuous realtime monitoring of polymerization reactions

#### Abstract

ACOMP (Automatic Continuous Online Monitoring of Polymerization Reactions) is a patented product from Fluence Analytics that continuously monitors and characterizes polymers during polymerization or post-polymerization processing steps of laboratory or industrial scale polymer reactions. Beyond allowing the user to monitor polymerizations in real time, ACOMP enables the user an ability to control specific desired polymer characteristics.

During the 1980s Professor Wayne Reed, Fluence Analytics' CSO, began laying the research foundation with his work in light scattering to characterize macromolecules while working at the Physics Department at Tulane University. His eventual goal was to analyze the evolution of these macromolecular processes in real time, and in the process, he developed a methodology capable of studying many polymerization techniques such as free radical, controlled radical, solution, bulk and emulsion during synthesis. Over the following 20 plus years, a multitude of government and industrially funded research projects and collaborations have led to over 90 invited lectures and over 40 peer review publications. Additionally, there have been 10 issued and 25 pending patents around the space of polymerization monitoring and control.

Today, ACOMP, as offered by Fluence Analytics, is a completely automated solution which allows for tighter process control, resulting in increased polymer yield, increased polymer quality and consistency, increased worker safety, reduced cycle time, reduced VOCs and reduced energy and raw material usage.

This technical note will explain how ACOMP functions as well as some of the measurable benefits gained from utilizing the ACOMP technique.



Figure 1: Industrial ACOMP



1

# **ACOMP** Principle

Rather than placing probes directly into the polymerization reactor or slip streams, ACOMP continuously extracts a small stream of reactor contents during the process (typically 0.10-1.0 cm<sup>3</sup>/minute). Next, the extracted stream is immediately diluted and homogenized, through proprietary sample handling technology, with an appropriate solvent halting further polymerization. The conditioned stream of diluted polymer passes through a detector train which enables realtime characterization of many polymerization processes. The detectors, of which any or all may be used, are comprised of Ultra-Violet/Visible Absorption (UV/Vis), Differential Refractometer (DRI), dilute solution single capillary viscometer (Visc) and multi-angle light scattering (MALS), and others. The appropriate configuration of detectors is selected according to the desired polymer properties and chemistry to be monitored. The subsequent data are then used to calculate the Weight Average Molecular Weight (Mw), Intrinsic Viscosity (IV), monomer conversion kinetics, residual monomers, comonomer composition and drift, and other polymer properties traditionally measured through discrete, off-line analysis.

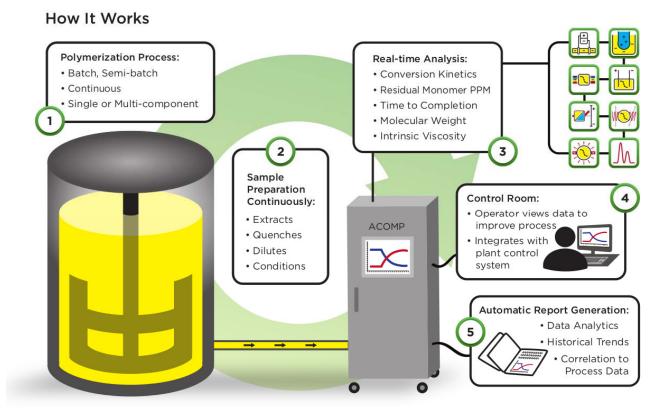


Figure 2: ACOMP process



# **Theoretical Framework**

Fluence Analytics can tailor ACOMP to pair with many detectors through a selection of modules. As noted, Fluence Analytics offers ACOMP with UV/Vis, DRI, Visc. and MALS. Other detectors that can be utilized in ACOMP include conductivity, pH, Oxygen sensing and near-infrared.

The UV/Vis detector uses multiple wavelengths whose selection is based on the evaluation of the absorption spectra of the monomer, polymer and solvent. By monitoring the absorption of UV light for specific wavelengths throughout the polymerization process, ACOMP can determine the monomer concentration and polymer conversion as well as actively predict, the total amount of residual monomer concentration.

The DRI detector is used to determine the total concentration of monomer and polymer in the sample stream.


Intrinsic Viscosity as measured by the Viscometer is a measure of both the hydrodynamic volume and molar mass of a polymer and can be correlated to the polymer's architecture and rheological properties. The intrinsic viscosity information can then be used to determine correlations related to the degree of polymer branching, crosslinking and shear thinning.

The Fluence Analytics built MALS detector is specifically designed for ACOMP use and utilizes multiangle Zimm extrapolation in conjunction with polymer concentration to characterize the weight average molecular weight of the polymer sample throughout the polymerization reaction. This information can be used to directly show the evolving molecular weight of the product as a function of time or polymer concentration.

# **Raw UV, Viscometer and Light Scattering Signals**

Figure 3 displays typical raw ACOMP signals versus time for the free radical polymerization of acrylamide. A monomer baseline is collected for all detectors prior to initiating the polymerization which is indicative of the known reactor conditions before any signal changes due to the effects of polymerization. The signal in red represents the UV Absorption versus time and is a direct measure of the monomer concentration throughout the polymerization. Subsequently, polymer concentration is calculated from monomer concentration through a simple mass balance. Also shown in figure 3 are 90-degree light scattering signature from the MALS detector and the raw voltage signal from the Viscosity detector. These detectors are molar mass sensitive and as result, increase throughout the polymerization as the size and concentration of the polymer increase. They are used to calculate the weight average molecular weight and intrinsic viscosity, respectively.

Figure 3: Typical online Raw UV/VIS, Light Scattering and Viscosity Signals for a batch reaction monitored by ACOMP

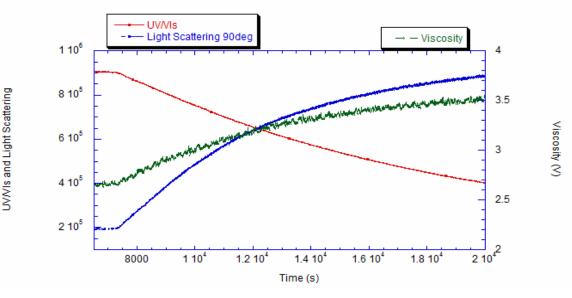



#### **Monomer and Polymer Concentration**

Figure 4 shows the calculated evolution of monomer and polymer concentrations throughout the polymerization. The consumption of monomer in a free radical polymerization follows first-order kinetics as the monomer double bonds form polymer bonds. Fluence Analytics has utilized this continuous data to develop a monomer prediction algorithm that can predict residual monomer concentration beyond the direct sensitivity of the UV/Vis detector. The results of this algorithm have been thoroughly cross-checked and confirmed via High Performance Liquid Chromatography (HPLC).

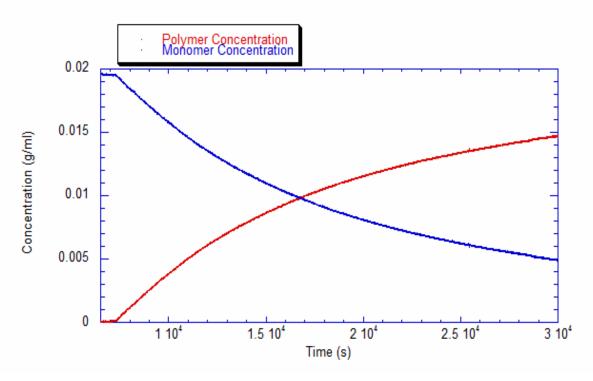


Figure 4: First order consumption of acrylamide and generation of polyacrylamide

#### **Intrinsic Viscosity and Molecular Weight**

As the polymer concentration is determined in real time, the weight average Molecular Weight M<sub>w</sub> and Intrinsic Viscosity are correspondingly calculated. Figure 5 shows the IV and M<sub>w</sub> for a free radical polymerization, where the kinetic chain length is proportional to the monomer concentration. Since monomer concentration decreases throughout the reaction, longer polymer chains are produced at the beginning of the reaction, and chain lengths diminish as the reaction proceeds. The ability to monitor the cumulative weight average molecular weight and polymer concentration also allows for the calculation of Instantaneous Weight Average Molecular Weight, M<sub>w,inst</sub>. The M<sub>w,inst</sub> measures the molecular weight of chains being produced at any instant, whereas the cumulative molecular weight applies to all the polymer chains produced up to any instant. Using these data, Fluence Analytics has developed a method for building "chromatography-free molecular weight distributions" by building the final distribution from the evolution of the instantaneous molecular weight throughout the reaction.



4

#### **Fluence Analytics**

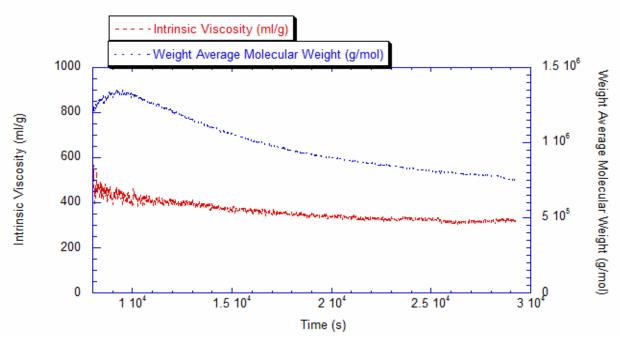


Figure 5: Molecular Weight and Intrinsic Viscosity

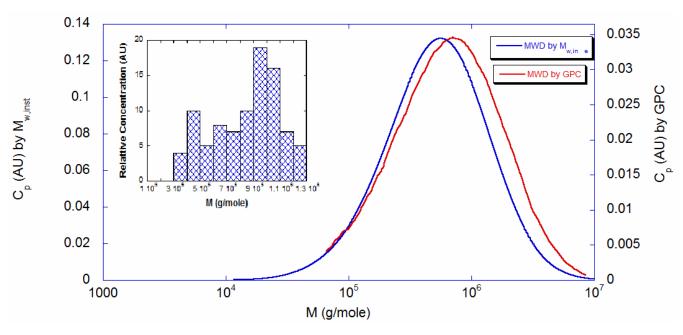


Figure 6: Molecular weight distribution by ACOMP M<sub>W,inst</sub> and GPC



# **ACOMP** Advantage

A major feature of ACOMP is the fully integrated and automated platform providing continuous, realtime analysis of polymer properties. This unique platform leverages the latest automation, sensor and detector technologies connected by a sophisticated hardware and software automation package. Data handling and advanced computational analysis packages have been developed to convert the continuous stream of ACOMP data into the critical information desired by the user. ACOMP's advantage is derived from direct polymer characterization continually taking place in real time, therefore, offering the capability of live feedback control.

Additionally, ACOMP has the ability to characterize the polymer without using size exclusion chromatography (SEC). While the SEC process is a valuable, post-production method of characterization, it inherently creates a delay from sample collection to process critical result and only provides discrete data at specified intervals. This leaves gaps in what is usually valuable reaction kinetic and material property evolution data during synthesis. These challenges, along with maintenance requirements, can make it difficult to adopt SEC directly to a rapidly changing industrial reaction process. ACOMP permits the continuous, realtime characterization of the polymer's absolute molecular weight, conversion of monomer, and intrinsic viscosity allowing for better monitoring and predictability of product characteristics. This translates to further control of the outcome of the polymerization process.

Increased Worker Safety	Reduced Cycle Time
Increased Product Quality	Reduced VOCs and Pollution
Increased Polymer Yield	Reduced Feedstock Usage
Increased Process Consistency	Reduced Energy Consumption
	Increased Product Quality Increased Polymer Yield

Figure 7: Measurable benefits of ACOMP usage

# Lab and Pilot ACOMP (IpACOMP)

The IpACOMP is available for laboratory and pilot scale polymerizations independently or customized to attach to, and monitor, pre-existing laboratory and/or pilot reactor systems. The ACOMP can also integrate with most standard control interfaces used by the laboratory or pilot reactors.

# Industrial ACOMP (iACOMP)

The Industrial ACOMP is a robust monitoring system designed for production-scale polymer reactors and can be tailored for numerous applications from a standardized set of modules.

Fluence Analytics' first Industrial ACOMP has been in service monitoring a batch specialty polymer process for more than two years with an overall availability greater than 97%. This availability is achieved by the use of more than 100 self-monitoring sensors coupled with automated corrective action procedures that engage as a result of any unplanned operating anomalies such as changes in system pressures. The Industrial ACOMP is designed in a NEMA 4X enclosure and can be rated for up to a Class I Division II plant environment. Class I Division I (explosion proof) environments can be considered with enclosure customization.

The ACOMP data are available to operators, plant managers, QA/QC personnel, process chemists and managers through the plant Distributed Control System (DCS).

Furthermore, iACOMP is capable of sending automated e-mail reports on each production batch or a specified period of time. These data can provide insights into Golden Batch information, the series of



#### Fluence Analytics

conditions that will yield on-spec product in the shortest amount of time, eliminate failed reactions, reduce grade change over time and off-spec product, and reduce batch cycle time and emissions.

### Conclusion

Fluence Analytics' ACOMP is the ONLY fully automatic, smart manufacturing system which enables the continuous, realtime monitoring and characterization of polymers during synthesis. Current characterization methods rely on off-line, single point characterization techniques which create significant information gaps throughout a reaction process. These industry standard techniques work very well for analyzing end products, but do not yield critical kinetic and property evolution information that is only available from ACOMP.

With this ACOMP enabled direct reaction monitoring capability, the next steps are being taken to include process optimization and fully automated reaction control. These continuous, realtime data can be fully integrated with control systems and other data analysis tools, thus creating another opportunity for long-term data mining and multi-site, coordinated product optimization. Finally, this level of control can enable tailored polymeric materials with customized properties and applications for individual customers.

#### For further reading about ACOMP:

Reed, W.F. "Automated Continuous Online Monitoring of Polymerization Reactions (ACOMP) and Related Techniques." Encyclopedia of Analytical Chemistry, 2013, 1-40, DOI: 10.1002/9780470027318.a9288.

Alb, A. M.; Reed, W. F. "Fundamental Measurements in Online Polymerization Reaction Monitoring and Control with a Focus on ACOMP", Macromolecular Reaction Engineering, 2010, 4(8), 470-485, DOI: 10.1002/mren.200900079

Fluence Analytics 1078 South Gavoso Street New Orleans, LA 70125 USA T: 1 - 504 - 777 - 2805 F: 1 - 504 - 777 - 2818 Technical Note: ACOMP 001 © 2017 by Fluence Analytics All rights reserved.



Fluence Analytics' mission is to optimize efficiency and quality in the production of polymeric materials and biopharmaceuticals bv deploving innovative products leveraging continuous, realtime data streams.







LOUISIANA

ENTERTAINMENT





**Our Partners** 







NewOrleansBioInnovationCenter

